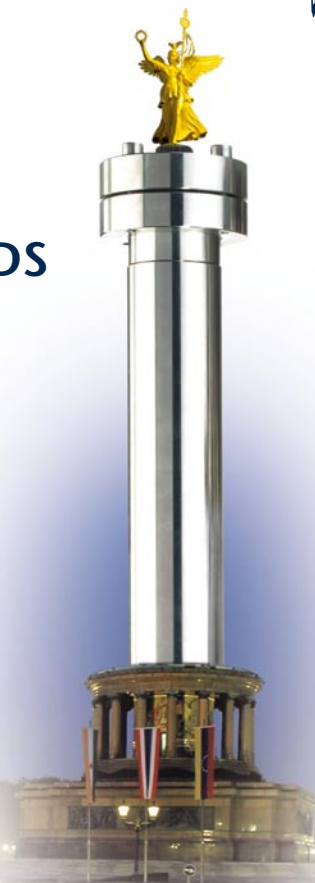
Euroline

ProntoSIL Hypersorb ODS

ProntoSIL Hypersorb ODS – the Alternative to Hypersil ODS

ProntoSIL Hypersorb ODS is our recommended alternative to Hypersil ODS. As with the original Hypersil ODS, our ProntoSIL Hypersorb ODS material is made for a broad range of applications. Like Hypersil ODS, ProntoSIL Hypersorb ODS is based on a first generation (type A) silica gel and has been developed to exhibit the same bonding density. Due to the slightly lower acidity of the basic silica gel, its silanophilic activity is a little bit lower than that of Hypersil ODS. The polarity of the ProntoSIL Hypersorb ODS surface, and hence also the selectivity of the material, is practically identical to that of Hypersil ODS. The same selectivity is also observed with ProntoSIL Hypersorb ODS although basic compounds do elute with better peak shape.

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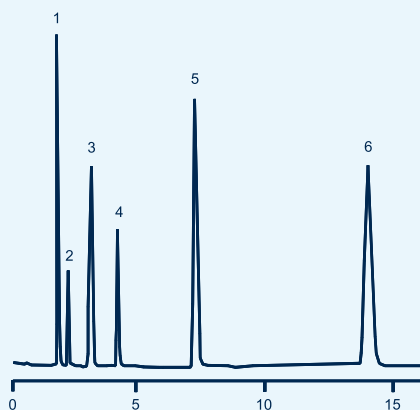
Hypersil ODS

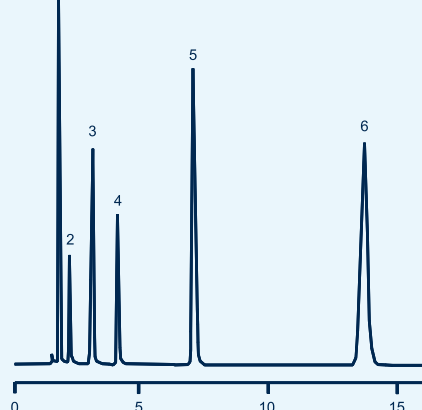
Test Mixture

1. benzamide
2. benzyl alcohol
3. acetophenone
4. methyl benzoate
5. benzophenone
6. biphenyl

Column: 150 x 4.6 mm
 Eluent: ACN/50 mM phosphate buffer (pH 3)
 Flow rate: 1.0 ml/min
 Temperature: 30 °C
 Detection: UV, 254 nm
 Inj. volume: 10 µl

ProntoSIL Hypersorb ODS

Benefits of replacement

- No revalidation of your HPLC method
- Same selectivity as with Hypersil ODS
- Comparable hydrophobicity
- Identical metal content
- Lower silanophilic activity
- Excellent value

Over 30 years of experience in the field of separation and extensive quality control ensure superior quality and the highest reproducibility of your analytical results. We design and manufacture HPLC column hardware ranging from 2 mm ID to 62 mm ID under strict quality control. We offer a wide range of column lengths from 5 mm up to 300 mm. An easily exchangeable integrated precolumn is available upon request.

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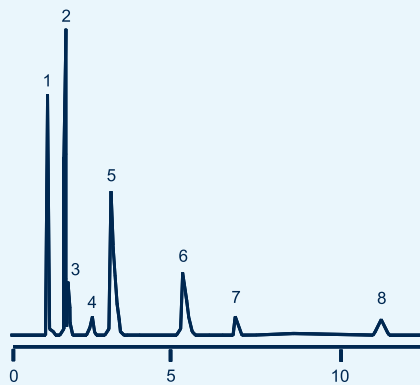
Quality Control

Every column that leaves our production has been individually checked and is shipped with a test certificate. We guarantee the quality of every HPLC column we deliver. Our performance standards with regard to peak symmetry and plate counts are extremely high to ensure consistent material and column quality.

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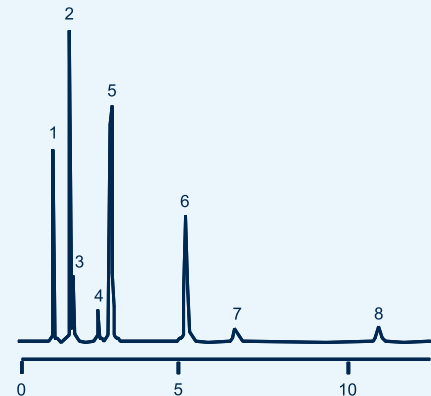
Column: 150 x 4.6 mm
 Eluent: MeOH/H₂O 49:51 (w/w)
 Flow rate: 1.0 ml/min
 Temperature: 40 °C
 Detection: UV, 254 nm
 Inj. volume: 5 µl

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Other column sizes are available upon request.
 Technical data are subject to change without notice.

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Visit www.knauer.net for details on complete HPLC systems, HPLC columns, and osmometers.

KNAUER · ASI · Advanced Scientific Instruments
 Wissenschaftliche Gerätebau
 Dr. Ing. Herbert Knauer GmbH
 Hegauer Weg 38 · D-14163 Berlin, Germany
 Telephone: +49 (0)30 / 80 97 27-0
 Telefax: +49 (0)30 / 8 01 50 10
 E-Mail: info@knauer.net
 Internet: www.knauer.net

Your local distributor: