Preparative HPLC Pump 1800

Reliable Power

KNAUER introduces the high performance **Preparative HPLC Pump 1800**.

It is the smallest preparative HPLC pump in its performance class with flow rates of up to 1,000 ml/min. It is based on latest engineering plus 30 years experience in developing HPLC pumps. All mechanical components are designed for continuous operation. Therefore the pump is also suitable for other precise dosing applications under high back pressures.

The impressive small housing of the **Preparative HPLC Pump 1800** gives you more degree of freedom for your preparative system setup with advantages like short fluid paths or small overall system size.

With integrated active pulsation dampening the **Preparative HPLC Pump 1800** delivers entirely the liquid phase without residual pulsation.

Four pump heads are available with flow rates of 100, 250, 500, or 1,000 ml/min respectively. Pump heads can easily be changed by loosening two capillaries and four screws. A cooling device to be mounted on the pump head allows operating even with liquid CO₂. A biocompatible version with titanium inlays is also available.

The preparative **Preparative HPLC Pump 1800** is characterized by long durability and low maintenance requirements. The pump head is service friendly designed so that it can easily be disassembled for validation or simple maintenance tasks like cleaning, changing of piston seals etc.

The **Preparative HPLC Pump 1800** demonstrates its flexibility by being upgradeable to as well high as low pressure gradient systems. It is controllable through ChromGate® or EuroChrom® software as well as from the integrated touch pad or through an open communication protocol.

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Technical Data and Ordering Information

Flow Properties

Delivery System Double Piston Pump

Max. Operating Pressure		Flow Rate Range
100 ml pump head	400 bar	0.1 – 99.9 ml/min
250 ml pump head	200 bar	0.1 – 249.9 ml/min
500 ml pump head	100 bar	0.1 – 499.9 ml/min
1,000 ml pump head	50 bar	0.1 – 999.9 ml/min

Flow Rate Precision

± 2% within 5–50% of flow range; Methanol/H₂O 80/20, v/v

Reproducibility RSD ≤ 0.1%

Pulsation Electronic compensation of residual pulsation

Gradients

High pressure gradient: Configurable master pump system with up to 3 slave pumps

Low pressure gradient: Control of up to 4 solenoid valves

Technical Parameters

Display 2x24 digits

System Protection Soft start, P_{min} and P_{max} adjustable

Control 2 RS-232 interfaces, 2 remote control strips, touchpad

Programming

max. 20 programs (1 configurable as emergency program) + 9 program links

GLP Function detailed report

Firmware Update downloadable by a service technician

Power Supply 115/230 V, 47–63 Hz, 450 W

Dimensions/Weight 220 x 185 x 340 mm (W x H x D) / 10.1 kg

Optional Accessories

Pump heads (with various inlays), cooling and heating device for pump heads, low pressure gradient valve blocks (ternary) up to 220 ml/min and (binary) up to 1,000 ml/min, inlet merging device, 8 channel pump head (8 x 0.1 – 50.0 ml/min, 200 bar)

Ordering Information

Order No.	Preparative HPLC Pump 1800
A40209-1	incl. 100 ml/min pump head with stainless steel inlays
A40201-1	incl. 250 ml/min pump head with stainless steel inlays
A40205-1	incl. 500 ml/min pump head with stainless steel inlays
A40202-1	incl. 1,000 ml/min pump head with stainless steel inlays

Order No.	Accessories
A4099-1	Low pressure gradient, ternary, stainless steel up to 220 ml/min
A4094-1	Low pressure gradient, binary, stainless steel up to 1,000 ml/min
A4036	8 channel pump head
A1117	Inlet merging device

Technical data are subject to change without notice.

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Dr. Ing. Herbert Knauer GmbH